


Vernier Go Wireless[®] Charging Station

Order Code GW-CRG

Charging Instructions

Connect one or more Go Wireless devices or supported Go Direct[™] sensors to the Go Wireless Charging Station for two hours.

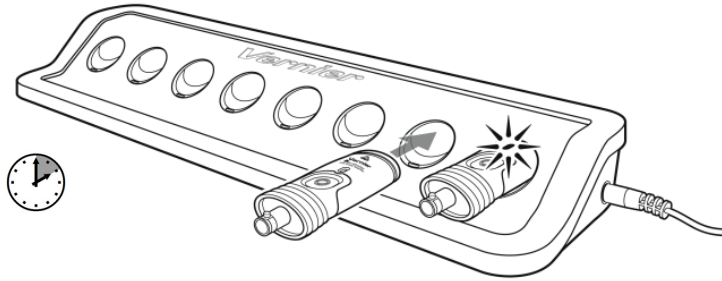

Blue LED Indicator on the Wireless Device

On – charging in progress

Off – charging complete

Precautions and Operational Tips

- You can charge devices for any amount of time, even for a few minutes between classes.
- Never insert metal objects into the charging slots.
- The Charging Station is not water resistant and should not be exposed to a wet environment.
- Do not insert a wet device into the Charging Station.

Disposal Instruction


When disposing of this electronic product, do not treat it as household waste. Its disposal is subject to regulations that vary by country and region. This item should be given to an applicable collection point for the recycling of electrical and electronic equipment. By ensuring that this product is disposed of correctly, you help prevent potential negative consequences on human health or on the environment. The recycling of materials will help to conserve natural resources. For more detailed information about recycling this product, contact your local city office or your disposal service.

The symbol, shown here, indicates that this product must not be disposed of in a standard waste container.



Warranty

Vernier warrants this product to be free from defects in materials and workmanship for a period of five years from the date of shipment to the customer. This warranty does not cover damage to the product caused by abuse or improper use.

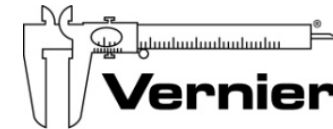

Measure. Analyze. Learn.[™]

Vernier Software & Technology

13979 S.W. Millikan Way • Beaverton, OR 97005-2886
Toll Free (888) 837-6437 • (503) 277-2299 • FAX (503) 277-2440
info@vernier.com • www.vernier.com

Rev 3/7/2017

Go Wireless is our registered trademark in the United States. Go Direct is our trademark in the United States.



Printed on recycled paper