

Go Direct[®] UV-VIS Spectrophotometer

(Order Code GDX-SPEC-UV)



The Go Direct UV-VIS Spectrophotometer is a small, lightweight ultraviolet and visible light spectrophotometer.

Note: Vernier products are designed for educational use. Our products are not designed nor are they recommended for any industrial, medical, or commercial process such as life support, patient diagnosis, control of a manufacturing process, or industrial testing of any kind.

What's Included

- Go Direct UV-VIS Spectrophotometer
- AC Power Supply
- Starna quartz cuvettes (2)
- Mini USB cable
- USB Power Adapter

Compatible Software and Interfaces

See www.vernier.com/manuals/gdx-spec-uv for a list of interfaces and software compatible with the Go Direct UV-VIS Spectrophotometer.

Quick Start: Vernier Spectral Analysis[®] and Bluetooth[®]

1. Connect the AC Power Supply to the spectrophotometer. This supplies power to the lamp. Turn the power switch to the ON position. Wait for the lamp indicator LED to remain green.
2. To power the Bluetooth radio, connect the spectrometer to the USB power adapter or to a powered USB hub. The Bluetooth[®] LED will blink. **Note:** Do not plug the USB into a computer USB port to run with Bluetooth wireless technology.
3. Launch Spectral Analysis, then click **Connect a Spectrometer**.
4. Select your sensor from the list. The sensor ID is located on the sensor label near the bar code. The Bluetooth LED on the sensor will now glow blue (no longer flashing).
5. Click **DONE**. You are now ready to collect data.

Using other Vernier data-collection apps or want to connect via USB?

Visit www.vernier.com/start-go-direct

Note: This sensor also works with LabQuest 2 and LabQuest 3; it does not work with the original LabQuest.

Powering the Spectrometer

Turning on the spectrometer	Connect to AC power. Turn the power switch on. Green LED indicator next to Power icon is solid when unit is ready to use. If indicator light turns red, disconnect all power, wait one full minute, and attempt again.
Turning off the spectrometer	Turn the power switch to the off position. The lamp will turn off and the fan will continue to run to cool the electronics for several more minutes. Wait until this is complete before disconnecting AC power.

Connecting the Spectrometer

See the following link for up-to-date connection information:

www.vernier.com/start/gdx-spec-uv

Connecting via Bluetooth

Ready to connect	Connect to AC power. Turn the power switch on. To power the Bluetooth radio, connect the spectrometer to the USB power adapter or to a powered USB hub. Blue LED next to Bluetooth icon flashes when sensor is ready to connect.
Connected	Blue LED next to Bluetooth icon is solid when sensor is connected via Bluetooth wireless technology.

Connecting via USB

Connected	LED next to Bluetooth icon is off. LED next to power is solid green. LED next to USB icon is solid green.
-----------	---

Using the Product with Spectral Analysis

Connect the sensor following the steps in the Quick Start section of this user manual.

Select the Type of Data to Measure

The three options for Experiment Types are

1. Measurement vs. Wavelength—collect a full spectrum.
2. Measurement vs. Concentration—conduct a Beer's law experiment.
3. Measurement vs. Time—collect time-based data for a kinetics experiment.

By default, Absorbance is expanded. You can also choose from % Transmittance, Emissions, or Advanced Full Spectrum. Use the Advanced Full Spectrum mode to switch among all the measurement types in the same session. For complete instructions, refer to the Spectral Analysis user manual at www.vernier.com/sa

Measurement vs. Wavelength (Full Spectrum)

1. Select Measurement vs. Wavelength.
2. If applicable, follow instructions for calibration. Align the cuvette so the clear side of the cuvette is facing the light source. **Note:** Calibration is not required for emissions.
3. You are now ready to collect data. Fill a cuvette about 3/4 full of a sample of the solution to be tested. Place the sample in the Spectrophotometer and click or tap Collect. Click or tap Stop to end data collection. The spectrum is automatically stored.
4. Save or export your data from the File menu.

Measurement vs. Concentration (Beer's law)

1. Select Measurement vs. Concentration.
2. If applicable, follow instructions for calibration. Align the cuvette so the clear side of the cuvette is facing the light source. **Note:** Calibration is not required for emissions.
3. Follow the instructions in the Choose a Wavelength dialog box. Select Done.
4. Click or tap Collect. Your first sample should still be in the Spectrophotometer. After the reading stabilizes, click or tap Keep. Enter the concentration of the sample and click or tap Keep Point.
5. Place your second sample in the cuvette slot. After the reading stabilizes, click or tap Keep. Enter the concentration of the sample and click or tap Keep Point.
6. Repeat the previous step for the remaining samples. When finished, click or tap Stop to end data collection. The data are automatically stored.
7. To see the best fit line equation for the standard solutions, click or tap Graph Tools, select Apply Curve Fit, and choose Linear. Click or tap Apply.
8. If doing Beer's law to determine the concentration of an unknown, place the unknown sample in the cuvette holder. Click or tap Graph Tools and enable Interpolate. Click or tap along the line until you find the concentration value that matches your unknown's measurement.
9. Save or export your data from the File menu.

Measurement vs. Time (Kinetics)

1. Select Measurement vs. Time.
2. If applicable, follow instructions for calibration. Align the cuvette so the clear side of the cuvette is facing the light source. **Note:** Calibration is not required for emissions.
3. Follow the instructions in the Choose a Wavelength dialog box. Select Done.
4. The default data-collection settings collect measurements every two seconds until the user manually stops data collection.
5. Mix the reactants. Transfer ~2 mL of the reaction mixture to a cuvette and place the cuvette in the spectrometer. Click or tap Collect.
6. When finished, click or tap Stop.
7. To fit a function to the data, click or tap Graph Tools, select Apply Curve Fit, and choose the appropriate curve fit. Click or tap Apply.
8. To add a calculated column to the data set, click or tap Column Options in the measurement header on the data table. Select Add Calculated Column. Modify

the name, units, and displayed precision accordingly. Select Insert Expression and select the appropriate equation. Modify the parameters and column options, if necessary. Click or tap Apply. The calculated column is automatically displayed on the graph.

9. Save or export your data from the File menu. To modify the data that are displayed on the graph, click or tap on the appropriate axis label.

Change the Settings in Spectral Analysis

1. Click or tap the gear to show the Spectrometer Settings dialog.
2. There are three parameters listed in the dialog box:
 - **Integration Time:** This is similar to the shutter speed of a camera. Spectral Analysis automatically selects the proper sample time during calibration in Absorbance and % Transmittance modes. If using Emissions or Advanced Experiment modes, you can modify this value to increase the amount of light hitting the detector and the signal spectral output.
 - **Wavelength Smoothing:** This is the number of adjacent readings on either side of a given value that is used to calculate an average value.
 - **Temporal Averaging:** This is the number of readings taken at a given wavelength to calculate an average reading.
3. Select the Calibrate button to recalibrate your spectrometer at any time.

Measure an Emission Spectrum with Spectral Analysis

You may use your Spectrophotometer to measure the visible emission spectrum of a light source such as an LED or a gas discharge tube. To do so, you will need to purchase the Vernier Optical Fiber (order code: VSP-FIBER).

Measure Intensity of Light Emissions

Insert the Vernier Optical Fiber into the spectrometer.

1. Connect the spectrometer following the steps in the Quick Start section of this user manual.
2. Launch Spectral Analysis.
3. Select the appropriate Emissions experiment from the listed options and follow the prompts in the app. Intensity is a relative measure with a range of 0–1. **Note:** The Spectrophotometer is not calibrated for measuring intensity.
4. Aim the tip of the optical fiber at a light source. Start data collection. Click or tap the Stop button to end data collection.

If the spectrum maxes out (flat and wide peaks at a value of 1), increase the distance between the light source and the tip of the optical fiber cable or reduce the integration time (see the Change the Settings in Spectral Analysis section).

To adjust the integration time, click or tap the gear. Set the integration time to a suitable value.

Using the Spectrophotometer with LabQuest

Select the Type of Data (or Units) You Want to Measure

There are three general types of data collection that measure absorbance or transmittance—absorbance (or %T) vs. wavelength, which produces a spectrum,

absorbance (or %T) *vs.* concentration for Beer's law experiments, and absorbance (or %T) *vs.* time for kinetics experiments.

The default data type is absorbance. If you want to measure the absorbance of a solution, proceed directly to the Calibrate section.

If you want to measure %T or intensity, do the following:

1. From the Sensors menu, choose Change Units ► USB: Spectrophotometer.
2. Select the unit or data type you wish to measure.

Calibrate the Spectrophotometer (Not Required if Measuring Intensity)

1. Choose Calibrate ► USB: Spectrophotometer from the Sensors menu.
Note: For best results, allow the Spectrophotometer to warm up for a minimum of five minutes.
2. Fill a cuvette about 3/4 full with distilled water (or the solvent being used in the experiment) to serve as the blank. After the Spectrophotometer has warmed up, place the blank cuvette in the Spectrophotometer. Align the cuvette so a clear side of the cuvette is facing the light source.
3. Follow the instructions in the dialog box to complete the calibration, and then tap OK.

Collect Data with LabQuest

Measurement vs. Wavelength (Generate a Spectrum)

1. Fill a cuvette about 3/4 full of the solution to be tested and place it in the Spectrophotometer.
2. Start data collection by tapping on the Start button in the lower left corner of the screen. Tap the Stop button to end data collection.
3. Select wavelength. You can tap on the graph to select a wavelength. Another way to change the wavelength is to navigate to the Meter screen, tap on the meter, and select Change Wavelength. Enter the wavelength of your choice and select OK. If the wavelength you type in is not measured by the unit, LabQuest will automatically choose the wavelength closest to your choice.
4. To store the spectrum data, tap on the file cabinet icon in the upper right of your screen.

Measurement vs. Concentration (Beer's Law Studies)

1. Generate a spectrum as described above. On the Meter screen, tap Mode. Change the mode to Events with Entry.
2. Enter the Name (e.g., Concentration) and Units (e.g., mol/L). Select OK.
3. A message will appear warning you to either save or discard the full spectrum run. Make your choice and proceed with data collection.
4. Place your first Beer's law standard solution in the Spectrophotometer. Start data collection. After the absorbance reading stabilizes, tap Keep. Enter the concentration of the solution and select OK.
5. Place your second standard sample in the Spectrophotometer. After the absorbance readings stabilize, tap Keep. Enter the concentration of the second sample and select OK.
6. Repeat Step 5 for the remaining standard samples. After you have tested the final standard, tap the Stop button to end data collection.

7. To calculate a best fit line equation for your standards, choose Curve Fit from the Analyze menu. Select Linear for the Fit Equation, and then select OK. The graph screen will appear again with the linear regression equation displayed.
8. Place a cuvette containing an unknown sample of solution in the Spectrophotometer. Tap the Meter tab and write down the displayed absorbance value. Tap the Graph tab and choose Interpolate from the Analyze menu. Trace the linear regression equation to determine the concentration of the unknown.

Measurement vs. Time (Kinetics)

1. Generate a spectrum as described above. On the Meter screen, tap Mode. Change the data-collection mode to Time Based.
2. You can change the rate, interval, and/or duration of time of data collection, if desired. Select OK when you are ready to proceed.
3. A message will appear warning you to either save or discard the full spectrum run. Make your choice and proceed with data collection.
4. Mix the reactants, transfer ~2 mL of the reaction mixture to a cuvette and place the cuvette in the Spectrophotometer. Start data collection. You may tap the Stop button to end data collection early.
5. To calculate a function for your data, choose Curve Fit from the Analyze menu. Select the Fit Equation, and then select OK. The graph screen will appear again.

Measure an Emission Spectrum with LabQuest

You may use your Spectrophotometer to measure the visible emission spectrum of a light source such as an LED or a gas discharge tube. To do so, you will need to purchase an optical fiber assembly (order code: VSP-FIBER).

Measure Intensity of Light Emissions

1. Insert the Spectrophotometer Optical Fiber into the UV-VIS Spectrophotometer.
2. Aim the tip of the optical fiber at a light source. Start data collection. Tap the Stop button to end data collection. **Note:** The Spectrophotometer is not calibrated for measuring intensity.

If the spectrum maxes out (flat and wide peaks at a value of 1), increase the distance between the light source and the tip of the optical fiber cable or reduce the sample time (see Change the Settings in LabQuest below).

To increase the sample time, or if data collection is unusually slow, choose Set Up Sensors ► Spectrophotometer: 1 from the Experiment menu. Set the Sample Time (begin with 75 ms, with subsequent reductions by 20 ms) to a suitable value and decrease the Samples to Average to 1.

Change the Settings in LabQuest

The Data Collection Screen in LabQuest lists all the settings for the device. To display this box choose Sensors ► Data Collection from the Meter screen.

For most experiments, the default settings work well.

There are five parameters listed in the dialog box.

- **Sample Time:** This is similar to the shutter speed of a camera. LabQuest automatically selects the proper sample time during calibration. **Note:** For emission studies, you may need to change the sample time manually.
- **Wavelength Smoothing:** This is the number of adjacent readings on either side of a given value that is used to calculate an average value. **Note:** Be careful adjusting this parameter as it may shift your wavelength values slightly.
- **Samples to Average:** This is the number of readings taken at a given wavelength to calculate an average reading.
- **Wavelength Range:** The range is determined by the type of Spectrophotometer in use.

Videos

View videos related to this product at www.vernier.com/gdx-spec-uv

Specifications

Detection modes	Absorbance, % Transmittance, Raw Lamp Output, and Intensity (fiber required)
Dimensions	18.5 cm × 17 cm × 7 cm
Lamp power supply	AC adapter (included)
Power consumption	3 A start-up, 500 mA continuous
Bluetooth power supply	USB (USB adapter included)
Absorbance light source	Deuterium (UV) and Incandescent (VIS)
Detector	Linear CCD
Wavelength range	220 nm–850 nm
Wavelength reporting interval	~1 nm
Optical resolution	3.0 nm (as determined with 486 nm hydrogen emission spectral line FWHM)
Wavelength accuracy	± 2.0 nm (as determined with holmium oxide NIST standard)
Absorbance photometric accuracy	± 5.0% (as determined with potassium dichromate NIST standards)
Typical scan time	~2 s
Sample format	10 mm × 10 mm cuvette (UV quartz cuvette included)
Beam height (z dimension)	8.5 mm

Safety

- Always wear UV blocking safety glasses or goggles when using this device. The light source emits ultraviolet radiation and can cause eye damage.

- When this instrument is in data-collection mode labeled Intensity, the light source will be blocked or turned off. Continue to use proper safety precautions.
- This instrument contains an internal high-voltage power source. Position the instrument where the ON/OFF switch can be easily reached. Do not move the instrument when it is running. Do not attempt to open or remove the case at any time.
- Do not remove or modify any of the installed safety components of this device. Doing so will create an unsafe operating condition and will void the product warranty.
- No user-serviceable parts are in this device. Do not attempt to open or modify this device. Contact Vernier for all repairs and service including lamp replacement.
- Handle the device with care. This instrument can be damaged if it is dropped.
- Do not use this instrument if it is damaged in any way. Contact Vernier Technical Support for troubleshooting and technical assistance.
- Do not use this instrument for clinical or diagnostic procedures.

Troubleshooting

1. For best results, the absorbance values for all samples must be between 0.1 and 1.0 absorbance units at all wavelengths.
2. Make sure you are using UV-VIS compatible cuvettes if you choose not to use the enclosed quartz cuvettes.

Find troubleshooting tips here: www.vernier.com/til/3044

Repair Information

- The UV light source in the Go Direct UV-VIS Spectrophotometer is a deuterium lamp. The lifetime of this source is rated at approximately 1000 hours. Lifetime is defined as the amount of time it takes the lamp to decrease to 50% of its original output. Deuterium lamp life is also inversely proportional to the number of ignitions, so it is best to limit the number of times you ignite the bulb if you wish to prolong its life.
- The deuterium lamp is covered by a one-year warranty or 1000 hours, whichever comes first.
- Contact Vernier for all repairs and service including lamp replacement
- No user-serviceable parts are in this device. Do not attempt to open the case of the device. Do not attempt to change or repair the lamp. Doing so will create an unsafe operating condition and will void the product warranty.

If you have watched the related product video(s), followed the troubleshooting steps, and are still having trouble with your Go Direct UV-VIS Spectrophotometer, contact Vernier Technical Support at support@vernier.com or call 888-837-6437. Support specialists will work with you to determine if the unit needs to be sent in for repair. At that time, a Return Merchandise Authorization (RMA) number will be issued and instructions will be communicated on how to return the unit for repair.

Accessories/ Replacements

Item

UV Quartz Cuvettes (2)

Spectrophotometer Optical Fiber

Mini USB Cable

Replacement Power Supply for Ultraviolet Range (UV-VIS) Spectrophotometers

Order Code

CUV-QUARTZ

VSP-FIBER

CB-USB-MINI

VSP-UV-PS

Warranty

Warranty information for this product can be found on the Support tab at www.vernier.com/gdx-spec-uv

General warranty information can be found at www.vernier.com/warranty

Federal Communication Commission Interference Statement

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

Reorient or relocate the receiving antenna.

Increase the separation between the equipment and receiver.

Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.

Consult the dealer or an experienced radio/TV technician for help.

FCC Caution

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:

(1) this device may not cause harmful interference and

(2) this device must accept any interference received, including interference that may cause undesired operation

RF Exposure Warning

The equipment complies with RF exposure limits set forth for an uncontrolled environment. The antenna(s) used for this transmitter must not be co-located or operating in conjunction with any other antenna or transmitter. You are cautioned that changes or modifications not expressly approved by the party responsible for compliance could void your authority to operate the equipment.

IC Statement

This device complies with Industry Canada license-exempt RSS standard(s). Operation is subject to the following two conditions:

(1) this device may not cause interference, and

(2) this device must accept any interference, including interference that may cause undesired operation of the device.

Industry Canada - Class B This digital apparatus does not exceed the Class B limits for radio noise emissions from digital apparatus as set out in the interference-causing equipment standard entitled "Digital Apparatus," ICES-003 of Industry Canada. Operation is subject to the following two conditions: (1) this device may not cause interference, and

(2) this device must accept any interference, including interference that may cause undesired operation of the device.

To reduce potential radio interference to other users, the antenna type and its gain should be so chosen that the equivalent isotropically radiated power (e.i.r.p.) is not more than that permitted for successful communication.

RF exposure warning: The equipment complies with RF exposure limits set forth for an uncontrolled environment. The antenna(s) used for this transmitter must not be co-located or operating in conjunction with any other antenna or transmitter.

Le présent appareil est conforme aux CNR d'Industrie Canada applicables aux appareils radio exempts de licence. L'exploitation est autorisée aux deux conditions suivantes :

(1) l'appareil ne doit pas produire de brouillage, et

(2) l'appareil doit accepter toute interférence radioélectrique, même si cela résulte à un brouillage susceptible d'en compromettre le fonctionnement.

Cet appareil numérique respecte les limites de bruits radioélectriques applicables aux appareils numériques de Classe B prescrites dans la norme sur le matériel interférent-brouilleur: "Appareils Numériques," NMB-003 édictée par Industrie Canada. L'utilisation est soumise aux deux conditions suivantes:

(1) cet appareil ne peut causer d'interférences, et

(2) cet appareil doit accepter toutes interférences, y comprises celles susceptibles de provoquer un dysfonctionnement du dispositif.

Afin de réduire les interférences radio potentielles pour les autres utilisateurs, le type d'antenne et son gain doivent être choisis de telle façon que l'équivalent de puissance isotrope émise (e.i.r.p.) n'est pas plus grand que celui permis pour une communication établie.

Avertissement d'exposition RF: L'équipement est conforme aux limites d'exposition aux RF établies pour un environnement non supervisé. L'antenne (s) utilisée pour ce transmetteur ne doit pas être jumelée ou fonctionner en conjonction avec toute autre antenne ou transmetteur.

Note: This product is a sensitive measurement device. For best results, use the cables that were provided. Keep the device away from electromagnetic noise sources, such as microwaves, monitors, electric motors, and appliances.



Vernier Science Education

13979 SW Millikan Way • Beaverton, OR 97005-2886

Toll Free (888) 837-6437 • (503) 277-2299 • Fax (503) 277-2440

info@vernier.com • www.vernier.com

Rev. 7/10/2024

Go Direct, Vernier Spectral Analysis, LabQuest, and other marks shown are our trademarks or registered trademarks in the United States.

All other marks not owned by us that appear herein are the property of their respective owners, who may or may not be affiliated with, connected to, or sponsored by us.

