

Go! Link[®]

(GO-LINK)




Go! Link is a single-channel USB interface used to connect many Vernier sensors directly to a Windows[®] or Macintosh[®] computer. Some Vernier sensors can also be used when Go! Link is connected to a Chromebook[™]. For a complete list of compatible sensors, visit www.vernier.com/go

Getting Started

Go! Link with a Computer

1. Install Logger Lite[®] or Logger Pro[®] software on your computer.
 - Logger Lite is free and can be downloaded at www.vernier.com/loggerlite
 - Logger Pro has many additional features and is available for purchase at www.vernier.com/lp
2. Connect Go! Link to the computer's USB port and the sensor to the Go! Link.
3. Start the software. The sensor will be recognized and in most cases will display the live reading. Note: The thermometer shown below is only displayed when using a temperature probe.
4. Click Collect to begin collecting data, if desired, or open an experiment.

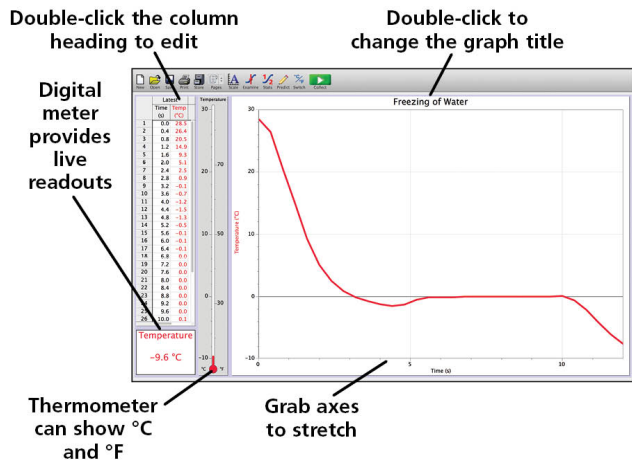

Go! Link with a Chromebook

1. Install Graphical Analysis[™] (version 1.2 or newer, available in the Chrome web store in late spring 2015).
2. Connect Go! Link to the Chromebook's USB port and the sensor to the Go! Link.
3. Launch Graphical Analysis. The sensor will be recognized and in most cases will display the live reading.
4. Start data collection, if desired.

NOTE: Vernier products are designed for educational use. Our products are not designed nor are they recommended for any industrial, medical, or commercial process such as life support, patient diagnosis, control of a manufacturing process, or industrial testing of any kind.

Specifications

USB specification	1.1
Maximum sampling rate	200 samples/second
Resolution	12 bit

Related Products

Go! Link Teacher Pack (order code: GL-TP)

Save money by purchasing a Teacher Pack of eight Go! Links. Includes free download of Logger Lite software. More information is available at www.vernier.com/go

Go! Temp[®] (order code: GO-TEMP)


Go!Temp is an easy-to-use temperature probe that connects directly to the USB port of a computer, LabQuest, or Chromebook. Includes free download of Logger Lite software. More information is available at www.vernier.com/go-temp

Go! Motion[®] (order code: GO-MOT)

Go! Motion is a motion detector that connects directly to the USB port of a computer, LabQuest, or Chromebook. Includes free download of Logger Lite software. More information is available at www.vernier.com/go-mot

Warranty

Vernier warrants this product to be free from defects in materials and workmanship for a period of five years from the date of shipment to the customer. This warranty does not cover damage to the product caused by abuse or improper use.

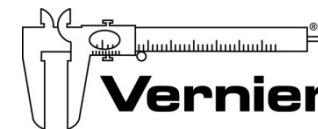

Measure. Analyze. Learn.[™]
Vernier Software & Technology

13979 S. W. Millikan Way • Beaverton, OR 97005-2886
Toll Free (888) 837-6437 • (503) 277-2299 • FAX (503) 277-2440
info@vernier.com • www.vernier.com

Rev. 1/12/15

Go! Link, Logger Pro, Logger Lite, Vernier LabQuest, Graphical Analysis, Go! Link, Go! Motion, and other marks shown are our trademarks or registered trademarks in the United States.

All other marks not owned by us that appear herein are the property of their respective owners, who may or may not be affiliated with, connected to, or sponsored by us.



Printed on recycled paper.

