

## Digital Adapter (Order Code DG-BTD)



This DG-BTD is used to adapt Vernier photogates, PASCO photogates, a PASCO Smart Pulley™, or the Radiation Monitor to LabQuest® 2, LabQuest, LabPro®, SensorDAQ®, or CBL 2™. Connect the quarter-inch stereo plug of the sensor cable to the adapter. Connect the white British Telecom plug of the adapter to one of the digital channels on the interface.

Please note that this adapter is not designed for use with the original CBL™. If you want to use photogates or Smart Pulleys with CBL, use the Photogate Adapter for CBL (order code CBL-P), or the Dual Photogate Adapter (order code CBL-2P).



Measure. Analyze. Learn.™

**Vernier Software & Technology**

13979 S.W. Millikan Way • Beaverton, OR 97005-2886

Toll Free (888) 837-6437 • (503) 277-2299 • FAX (503) 277-2440

info@vernier.com • www.vernier.com

Rev.9/19/2012