# Polar<sup>®</sup> Transmitter Module

## Order Code HR-TRANS

The Polar Transmitter Module is used with an Exercise Heart Rate Strap (HR-STRAP, not included) or Heart Rate Hand Grips (HR-GRIP, not included). Either combination is an excellent option for continuously monitoring heart rate before, during, and after exercise or while a person is stationary. Data are wirelessly transmitted to iPad<sup>®</sup>, LabQuest 2, or other mobile devices that are Bluetooth<sup>®</sup> Smart Ready. For information on how to use the Polar Transmitter Module with either of these products see the related user manuals at [www.vernier.com/gw-ehr](http://www.vernier.com/gw-ehr) and [www.vernier.com/gw-hr](http://www.vernier.com/gw-hr).

**NOTE:** Vernier products are designed for educational use. Our products are not designed nor are they recommended for any industrial, medical, or commercial process such as life support, patient diagnosis, control of a manufacturing process, or industrial testing of any kind.

## Specifications

### Polar Transmitter Module

Battery Type	CR 2025
Battery Lifetime	200 hrs
Operating Temperature	-10 to 50°C
Wireless Range	10 m or more unobstructed
Radios	Bluetooth and 5 kHz RF transmission

## Disposal Instruction

When disposing of this electronic product, do not treat it as household waste. Its disposal is subject to regulations that vary by country and region. This item should be given to an applicable collection point for the recycling of electrical and electronic equipment. By ensuring that this product is disposed of correctly, you help prevent potential negative consequences on human health or on the environment. The recycling of materials will help to conserve natural resources. For more detailed information about recycling this product, contact your local city office or your disposal service.

The symbol, shown here, indicates that this product must not be disposed of in a standard waste container.

## Warranty

Vernier warrants the Polar Transmitter Module to be free from defects in materials and workmanship for a period of two years from the date of shipment to the customer. This warranty does not cover the battery or damage to the product caused by abuse or improper use.

## Federal Communication Commission Interference Statement

This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to Part 15 of the FCC rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

### FCC Caution

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:

(1) this device may not cause harmful interference and

(2) this device must accept any interference received, including interference that may cause undesired operation

### RF Exposure Warning

The equipment complies with RF exposure limits set forth for an uncontrolled environment. The antenna(s) used for this transmitter must not be co-located or operating in conjunction with any other antenna or transmitter.

You are cautioned that changes or modifications not expressly approved by the party responsible for compliance could void your authority to operate the equipment.

### IC Statement

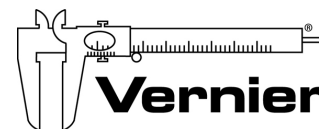
This device complies with Industry Canada license-exempt RSS standard(s). Operation is subject to the following two conditions: (1) this device may not cause interference, and (2) this device must accept any interference, including interference that may cause undesired operation of the device.

**Industry Canada - Class B** This digital apparatus does not exceed the Class B limits for radio noise emissions from digital apparatus as set out in the interference-causing equipment standard entitled "Digital Apparatus," ICES-003 of Industry Canada. Operation is subject to the following two conditions: (1) this device may not cause interference, and (2) this device must accept any interference, including interference that may cause undesired operation of the device. To reduce potential radio interference to other users, the antenna type and its gain should be so chosen that the equivalent isotropically radiated power (e.i.r.p.) is not more than that permitted for successful communication.

**RF exposure warning:** The equipment complies with RF exposure limits set forth for an uncontrolled environment. The antenna(s) used for this transmitter must not be co-located or operating in conjunction with any other antenna or transmitter.

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Measure. Analyze. Learn.<sup>™</sup>  
Vernier Software & Technology

13979 S. W. Millikan Way • Beaverton, OR 97005-2886

Toll Free (888) 837-6437 • (503) 277-2299 • FAX (503) 277-2440

[info@vernier.com](mailto:info@vernier.com) • [www.vernier.com](http://www.vernier.com)

Rev. 11/19/2020

The Polar Logo is a registered trademark of Polar Electro.

Printed on recycled paper.